


Comparator Kit Phosphate

PO_4 (0 - 80 mg/l)

1. Fill both cells with 10 ml. Place one cell in the left side of the comparator a blank.
2. Second cell: Add one tablet of *Phosphate HR* and crush with the stirring rod and mix gentle.
3. Insert the second cell with added regant tablet to the right side of the comparator.
4. Insert the Phosphate disc to the comparator and rotate it until the color match is obtained.
5. Now you can read the value.

Note!
If samples are turbid,
filter before testing.

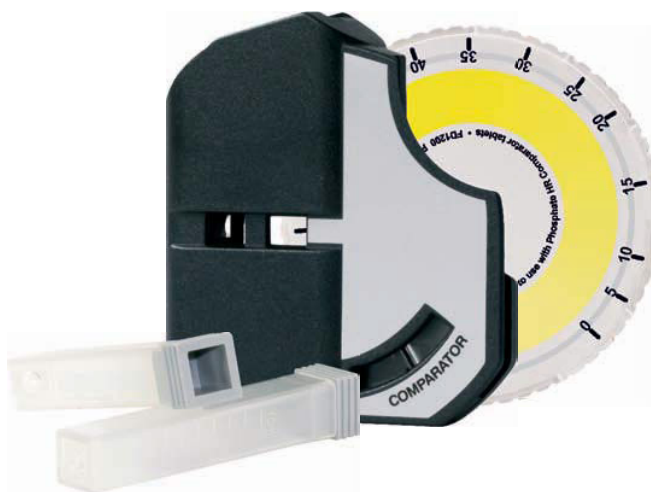

PCKFD1200

Kit contains:

Black plastic carrying case

Comparator PCP001

2 x 10 ml cell SPkv5

Stirring rod SPstr1

Cleaning brush SPclb1

Disc for Phosphate FD1200

50 tbl. Phosphates Comparator TbsRPP50



Attention!
Wear protective gloves and safety goggles when performing any tests using corrosive, harmful or irritant reagents. Do not ingest.